
MICROSCOPE & TELESCOPE KIT

Levenhuk LabZZ MT2

LabZZ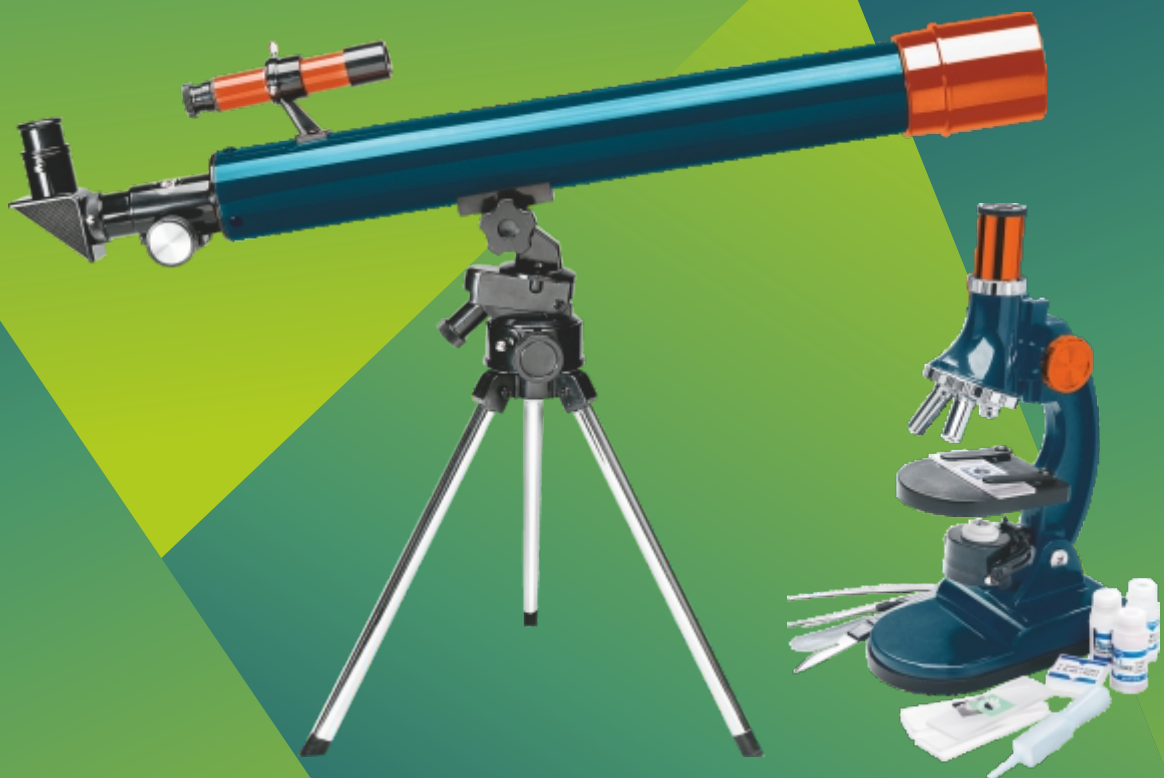

User Manual
Návod k použití
Bedienungsanleitung
Guía del usuario
Instrukcja obsługi
Инструкция по эксплуатации

levenhuk[®]
Zoom&Joy

Levenhuk LabZZ 75x/300x/900x Microscope

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Dear friend!

Levenhuk LabZZ MT2 Kit for children will open up the amazing world of exciting scientific discoveries! With the telescope in the kit you can study the Moon, planets and bright stars, unravel the mysteries of the endless Cosmos and observe the most distant objects. The microscope, on the other hand, will help you witness the unseen – things so small that you can't see them with the naked eye! The microscope in the kit will help you find out what things around us are made of and what kinds of secrets are hidden from us in every day life. Your experience with Levenhuk LabZZ MT2 Kit will be both informative and entertaining. Carefully read the entire manual before operating the telescope or microscope. Don't lose it; you might need it later.

Information for parents

Dear adults! Remember that you are responsible for your child at all times. Always ensure the child's safety when working with the microscope or telescope. Read this entire manual carefully, especially the handling and maintenance instructions. Even if your child is old enough and reads well, repeat all the rules aloud one more time before beginning to work with the instruments. Make sure that he or she clearly understands these instructions.

CAUTION! CHOKING HAZARD! These devices include small parts. The telescope and microscope are designed for children over 5 years of age and should only be used under adult supervision.

Care and maintenance (microscope and telescope)

- Never, under any circumstances, look directly at the Sun through this device without a special filter, or look at another bright source of light or at a laser, as this may cause PERMANENT RETINAL DAMAGE and may lead to BLINDNESS.
- You should ONLY observe the Sun with a professional solar aperture filter that completely covers the front of the instrument. The front end of the finderscope must be covered with aluminum foil or other non-transparent material to protect the internal parts of the telescope from damage. For safe solar observations we recommend using filters with an optical density (OD) of 5.0. Such filters transmit only 0.001% of the sun's light, blocking the remaining 99.999%. Most reliable solar filters are made of Baader AstroSolar film.
- Take necessary precautions when using the device with children or others who have not read or who do not fully understand these instructions.
- All parts of the telescope and microscope will arrive in one box. Be careful when unpacking it. We recommend keeping the original shipping containers. In the event that the telescope needs to be shipped to another location, having the proper shipping containers will help ensure that your telescope survives the journey intact. Make sure all the parts are present in the packaging. Be sure to check the box carefully, as some parts are small.
- Do not try to disassemble the device on your own for any reason. For repairs and cleaning of any kind, please contact your local specialized service center.
- Do not touch the optical surfaces with your fingers. To clean the device exterior, use only special cleaning wipes and special optics cleaning tools from Levenhuk. Do not use any corrosive or acetone-based fluids to clean the optics.
- Protect the device from sudden impact and excessive mechanical force. Do not apply excessive pressure when adjusting focus. Do not overtighten the locking screws.
- Store the device in a dry, cool place away from hazardous acids and other chemicals, away from heaters, open fire and other sources of high temperatures.
- If you are not using the device for extended periods of time, store the objective lenses and eyepieces separately from it.
- Be careful during your observations, always replace the dust cover after you are finished with observations to protect the device from dust and stains.
- Do not use the device for lengthy periods of time, or leave it unattended in direct sunlight. Keep the device away from water and high humidity.
- When using the microscope, try not to use it near flammable materials or substances (benzene, paper, cardboard, plastic, etc.), as the base may heat up during use, and might become a fire hazard.
- Always unplug the microscope from a power source before opening the base or changing the illumination lamp. Regardless of the lamp type (halogen or incandescent), give it some time to cool down before trying to change it, and always change it to a lamp of the same type.
- The chemicals might be harmful to human health. Use only those specified in the user manual. Make sure that your child knows how to use with them correctly.
- The workplace should be kept clean and tidy. Don't work with specimens in the kitchen near food you eat. The area should be well lit. You also should have fast and convenient access to clean water.
- Find a separate bucket for waste. All liquids should be poured down the drain, not into the sink.
- Dyes: methylene blue and methyl orange. Do not swallow. Keep out of the reach of children. If swallowed, seek medical advice immediately.
- Glue: usually organic material (for example, pitch). Do not swallow. Keep out of the reach of children. If swallowed, seek medical advice immediately.
- The kit includes everything needed for assembly, but if you need some additional tool you can use a basic set of standard instruments that almost everyone keeps at home.
- **Seek medical advice immediately if a small part or a battery is swallowed.**
- **Children should use the telescope under adult supervision only.**

Batteries safety instructions

- Always purchase the correct size and grade of battery most suitable for the intended use.
- Always replace the whole set of batteries at one time; taking care not to mix old and new ones, or batteries of different types.
- Clean the battery contacts and also those of the device prior to battery installation.
- Make sure the batteries are installed correctly with regard to polarity (+ and -).
- Remove batteries from equipment that is not to be used for an extended period of time.
- Remove used batteries promptly.
- Never attempt to recharge primary batteries as this may cause leakage, fire, or explosion.
- Never short-circuit batteries as this may lead to high temperatures, leakage, or explosion.
- Never heat batteries in order to revive them.

- Do not disassemble batteries.
- Remember to switch off devices after use.
- Keep batteries out of the reach of children, to avoid risk of ingestion, suffocation, or poisoning.
- Utilize used batteries as prescribed by your country laws.

Welcome to the microworld!

With Levenhuk LabZZ MT2 Microscope you will make lots of interesting discoveries and create exciting memories! Remember that following safety rules is absolutely necessary. In the kit you will find the microscope, as well as many useful accessories. Everything in place? Check it together with your parents. ④

Microscope's construction ①

1. **Eyepiece.** You observe the magnified image of the sample through the eyepiece. Eyepiece is a lens system that works the same way as a magnifying glass.
2. **Focusing knob.** Blurry images? Rotate the knob gently, adjusting the focus until all details become clear. Be careful - rotate the knob slowly so as not to damage the objective or observed slide.
3. **Monocular head.** This is a very important part of the microscope. The light travels right up the tube of the head directly into the eyepiece, where the resulting image is formed.
4. **Revolving nosepiece.** This is a device that holds three objectives with different magnifications. To change magnification, rotate the nosepiece until it clicks (the sound means that the chosen objective is in place).
5. **Objective.** This is needed for magnifying images. Levenhuk LabZZ MT2 Microscope comes with three objectives with different magnifications. They magnify images 75, 300 and 900 times compared to what you see with the naked eye.
6. **Stand.** This curved part of the microscope connects all other parts together. If you are feeling uncomfortable looking through the eyepiece, tilt or lift the stand until it's more comfortable. Change the inclination angle slowly, without abrupt movements.
7. **Stage.** This is exactly where you place the slide in order to observe it.
8. The stage has two **slide holders**, which, as is clear from their name, hold slides in place.
9. **The mirror** is located under the stage. It refracts the light from the lamp or the window illuminating the sample on the stage.
10. **Illumination lamp.** If illumination in your room is not enough, turn on the built-in illumination by turning over the mirror. ②
11. **Battery compartment.** The battery compartment is in the base of the microscope. What's it for? That's simple – it's where you

How to work with the microscope

1. First install batteries for illumination. Flip the microscope over and remove the cap from the battery compartment. Insert two AA batteries (those are not the thinnest penlight batteries, but those that are a bit thicker). Insert the batteries according to their polarity – make sure that the “+” sign on the battery matches “+” in the compartment (same goes for “-” signs). ③ Put the cap back on. Don't hesitate to ask adults for help if you are not sure.
2. Place the microscope on the flat surface near the source of bright light (window or a lamp). Turn the mirror towards the illumination source. You should see a bright light spot through the eyepiece.

CAUTION! Never direct the mirror towards the Sun, as it may ruin your eyesight and even cause blindness.

3. Now the microscope is ready for work.
4. Take a prepared microscope slide, place it on the stage and secure with holders.
5. Choose the objective by rotating the revolving nosepiece. The microscope comes with three objectives with different magnifications. The longer the objective, the higher magnification it produces. Most observations are better performed with lower magnification, so choose the shortest objective to start. You will hear a “click” when the objective is in place.
6. Slowly rotate the focusing knob so the objective is almost touching the slide. **BUT!** It shouldn't actually touch the slide as both the slide and the objective may be damaged if it does. That's why you need to do everything very slowly and carefully. When the objective is very close to the slide (but doesn't touch it!), look through the eyepiece and start to rotate the focusing knob in the opposite direction very slowly until you see a clear image.

Wonderful telescope

Telescope's construction ⑤

1. **Objective.** The most important part of the telescope. The objective is a lens system that gathers light from the distant object and forms its image. The objective should always be pointed in the direction of the object you want to observe.
2. **Optical tube.** Light rays collected by the objective travel down the optical tube to enter the eyepiece.
3. **Eyepiece.** You look at the Moon, planets and stars through the eyepiece. The eyepiece enlarges images of objects many times; that's why all objects look bigger than when observed with the naked eye.
4. **Sun shade.** The shade is a short tube that is put on the objective. It's needed to not let the light from street lamps, car lights, windows and other close-by light sources get in the objective and ruin the resulting image of the object you want to observe.
5. **Finderscope** – is a small telescope that is attached to the tube of the big telescope. The finderscope has small magnification and wide field of view (the latter means that you can see larger areas at once). The finderscope is made to help you find the desired object more quickly and easily. After you've found the object using the finderscope, you can observe it in detail through the telescope.
6. **Finderscope bracket** is required to attach the finderscope to the telescope tube.
7. **Focuser** – is a special device in which the eyepiece is inserted. The focuser is needed for bringing the eyepiece closer or farther from the objective. By doing so, you can achieve the maximum clarity of images.

8. If the image you see through the eyepiece isn't clear, rotate the **focusing knob** carefully.
9. **Diagonal mirror**. If you want to observe objects high above your head, use the diagonal mirror so you don't have to throw your head back. The diagonal mirror is inserted in the focuser. It deflects light rays so you can observe from a more comfortable position.
10. **Altazimuth mount** – is a special mechanism that allows you to move the telescope tube up-down and right-left.
11. **Azimuth lock knob** moves the tube horizontally.
12. **Altitude lock knob** moves the tube vertically.
13. **Slow-motion control**. The telescope is a very sensitive instrument; that's why for more precise pointing you need to use slow-motion control. This control allows you to move the tube very smoothly and without jumps.
14. **Tabletop tripod**. The telescope should be very stable so the images you see don't shake. This is why the tube is installed on a special support – the tripod.

How to assemble your telescope 6

It is very important to assemble the telescope correctly so it works the right way. Examine the telescope carefully. Try to remember the names of its parts, where each part goes and what it's needed for. It's better to assemble the telescope with adult help.

1. Open the box and put out all the parts. Double-check the box when it's empty – some small parts might still be in there.
2. Spread the tripod legs and install it on a flat surface so it is stable.
3. Find the mounting screw on the telescope (located underneath). Loosen it and set the telescope on the mount. Insert the screw in the holes on the telescope and the mount. Carefully tighten it. **Attention: Do not overtighten the screw as you may accidentally damage the screw thread.**
4. Loosen the finderscope mounting screws and slide it in the finderscope base. Carefully tighten the screws. The finderscope should be installed in a way so it points in the same direction as the telescope tube.

How to work with the telescope

Before starting observations it is very important to install eyepieces and all additional parts correctly.

Focuser: Turn the focusing knob toward the telescope until it will not go farther. Then start turning the knob towards yourself slowly until the image is clear. **Attention: Before focusing the view (adjusting its sharpness), make sure that eyepiece is inserted in the focuser. Otherwise, you won't see an image.**

18mm eyepiece: In the box with your telescope you will find two eyepieces. The lower the number on the eyepiece, the higher the magnification. It's better to start observing using an eyepiece with low magnification and wide field of view. The 18mm erecting eyepiece can be used for lunar and planetary observations and you can also observe a great number of distant terrestrial objects with it such as buildings, towers, antennas and others! **Attention: The 18mm erecting eyepiece should be inserted directly into the focuser, without the diagonal mirror.**

Diagonal mirror: Pull out the erecting eyepiece and insert the diagonal mirror into the focuser. 7

H12.5mm eyepiece: After the diagonal mirror is inserted into the focuser, insert the H12.5mm eyepiece in it. It provides higher magnification and is great for night-time observations.

How to start observing

**CAUTION! Never look at the Sun and the sky around it through a telescope!
It can ruin eyesight and even cause blindness!**

Before you start exploring the Cosmos, you should learn to operate the telescope during the day. Use the 18mm erecting eyepiece to observe different terrestrial objects – houses, trees, antennas on the rooftops and many others! This way you will learn to control the telescope and focus on desired objects. Besides, it's a lot of fun! **Attention: The telescope should be used in a place protected from the wind.**

When you get to observing the Moon, planets and stars at night, remember to choose locations away from street lamps, car lights and window lights. Try to observe on nights when the stars shine bright and evenly.

Point the telescope at the desired object, for example, the Moon. Looking through the finderscope, slowly move the tube until the object is in the center. Now look through the eyepiece and you will see the image of the object magnified many times!

When working with the telescope, be careful and patient. Don't push it and don't move it from its place. After some training you will learn to move the telescope tube without losing the object from the sight of the eyepiece.

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EN

Levenhuk International Lifetime Warranty

All Levenhuk telescopes, microscopes, binoculars and other optical products, except for accessories, carry a **lifetime warranty** against defects in materials and workmanship. Lifetime warranty is a guarantee on the lifetime of the product on the market. All Levenhuk accessories are warranted to be free of defects in materials and workmanship for **six months** from date of retail purchase. Levenhuk will repair or replace such product or part thereof which, upon inspection by Levenhuk, is found to be defective in materials or workmanship. As a condition to the obligation of Levenhuk to repair or replace such product, the product must be returned to Levenhuk together with proof of purchase satisfactory to Levenhuk.

This warranty does not cover consumable parts, such as bulbs (electrical, LED, halogen, energy-saving and other types of lamps), batteries (rechargeable and non-rechargeable), electrical consumables etc.

For further details, please visit our web site: <https://www.levenhuk.com/warranty>

If warranty problems arise, or if you need assistance in using your product, contact the local Levenhuk branch.

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Levenhuk prepared slides sets

The Levenhuk slides sets include thoroughly prepared specimens of various biomaterials as well as blank slides and cover glasses for you to make your own specimens. Prepared slides, blank slides and cover slips can be used with any microscope model.

Levenhuk Worldwide:

USA: www.levenhuk.com
Canada: ca.levenhuk.com
Czech Rep.: www.levenhuk.cz
Estonia: ee.levenhuk.com
Germany: de.levenhuk.com
Iceland: is.levenhuk.com

India: in.levenhuk.com
Latvia: lv.levenhuk.com
Lithuania: lt.levenhuk.com
Netherlands: nl.levenhuk.com
Poland: pl.levenhuk.com
Russia: www.levenhuk.ru

Slovakia: sk.levenhuk.com
Spain: es.levenhuk.com
Turkey: tr.levenhuk.com
Ukraine: www.levenhuk.ua
United Kingdom: uk.levenhuk.com
EU: eu.levenhuk.com

Levenhuk, Inc. 924-D East 124th Ave. Tampa, FL 33612 USA
Levenhuk®, LabZZ® are registered trademarks of Levenhuk, Inc.
© 2006–2017 Levenhuk, Inc. All rights reserved.
20161229

levenhuk[®]
Zoom&Joy[®]